



# Solutions Guide

DISINFECTANTS AND CLEANERS

TexQ™ | Bru-Clean TbC™ | TexCide™ | Hydrogen Peroxide | IPA | Ethanol

## Table of Contents

|                                          |    |
|------------------------------------------|----|
| TexQ Disinfectant .....                  | 2  |
| Bru-Clean TbC™ .....                     | 10 |
| TexCide™ .....                           | 16 |
| TexP Hydrogen Peroxide .....             | 22 |
| IPA .....                                | 26 |
| Denatured Ethanol .....                  | 30 |
| Crush Tube Swab .....                    | 32 |
| Appendix .....                           | 34 |
| Disinfectant Classification .....        | 35 |
| Disinfectant + Wiper Compatibility ..... | 36 |
| Pre-Wetted Wipers .....                  | 37 |
| ISO 5 Sample Cleaning Frequency .....    | 38 |



## About Texwipe

Texwipe's commitment to innovation, leadership and quality in cleanroom consumable products spans more than fifty years. We invest in technology to respond to our customers' evolving needs in contamination control.

### Texwipe's Core Values are:

- **Innovation** - Texwipe pioneers the latest technologies to provide innovation in contamination control products and processes.
- **Quality** - Texwipe product quality is maintained by the most advanced testing and quality control standards in the industry.
- **Technology Leadership** - Texwipe leads our industry in testing metrology, methods and processes to reduce contamination.

Throughout Texwipe's global operations, we support our customers with products designed to exceed the requirements for cleanroom consumable products. Our highly professional team will help you select and develop products for any critical environment application.



texwipe.com

# How to Choose

## Disinfectants/Cleaners Differentiation

|                                            | Disinfectants                                                                     |                                                                                   |                                                                                   | Cleaners                                                                            |                                                                                     |                                                                                     |
|--------------------------------------------|-----------------------------------------------------------------------------------|-----------------------------------------------------------------------------------|-----------------------------------------------------------------------------------|-------------------------------------------------------------------------------------|-------------------------------------------------------------------------------------|-------------------------------------------------------------------------------------|
|                                            |  |  |  |  |  |  |
|                                            | TexQ                                                                              | Bru-Clean                                                                         | TexCide                                                                           | TexP                                                                                | 70% IPA                                                                             | 70% Ethanol                                                                         |
| <b>Properties</b>                          |                                                                                   |                                                                                   |                                                                                   |                                                                                     |                                                                                     |                                                                                     |
| 0.2 micron filtered                        | ✓                                                                                 | n/a                                                                               | ✓                                                                                 | ✓                                                                                   | ✓                                                                                   | ✓                                                                                   |
| One-step cleaner/disinfectant              | ✓                                                                                 |                                                                                   | ✓                                                                                 |                                                                                     |                                                                                     |                                                                                     |
| Biodegradable                              |                                                                                   | ✓                                                                                 | ✓                                                                                 | ✓                                                                                   |                                                                                     |                                                                                     |
| EPA-registered disinfectant                | ✓                                                                                 | ✓                                                                                 | ✓                                                                                 |                                                                                     |                                                                                     |                                                                                     |
| No dyes or fragrances                      | ✓                                                                                 |                                                                                   | ✓                                                                                 |                                                                                     |                                                                                     |                                                                                     |
| No Volatile Organic Compounds (VOC)        | ✓*                                                                                |                                                                                   | 0.5% at use concentrate                                                           | ✓                                                                                   |                                                                                     |                                                                                     |
| Non-flammable                              | ✓                                                                                 | ✓                                                                                 | ✓                                                                                 | ✓                                                                                   |                                                                                     |                                                                                     |
| Shelf life, years                          | 2                                                                                 | 2                                                                                 | 1                                                                                 | 3                                                                                   | 2-3                                                                                 | 2                                                                                   |
| USP-compliant components (made with)       | n/a                                                                               | n/a                                                                               | n/a                                                                               | ✓                                                                                   | ✓                                                                                   | ✓                                                                                   |
| <b>Applications/Agents</b>                 |                                                                                   |                                                                                   |                                                                                   |                                                                                     |                                                                                     |                                                                                     |
| Cleaning                                   | ✓                                                                                 |                                                                                   | ✓                                                                                 | ✓                                                                                   | ✓                                                                                   | ✓                                                                                   |
| Disinfectant rotation program use          | ✓                                                                                 | ✓                                                                                 | ✓                                                                                 | ✓                                                                                   | ✓                                                                                   | ✓                                                                                   |
| Gamma-irradiated available                 | ✓                                                                                 |                                                                                   |                                                                                   |                                                                                     | ✓                                                                                   | ✓                                                                                   |
| Needs rinse                                | ✓                                                                                 | ✓                                                                                 |                                                                                   |                                                                                     |                                                                                     |                                                                                     |
| Residue removal agent                      |                                                                                   |                                                                                   |                                                                                   | ✓                                                                                   | ✓                                                                                   | ✓                                                                                   |
| <b>Testing</b>                             |                                                                                   |                                                                                   |                                                                                   |                                                                                     |                                                                                     |                                                                                     |
| Endotoxin tested                           | ✓                                                                                 |                                                                                   |                                                                                   |                                                                                     | ✓                                                                                   | ✓                                                                                   |
| Sterile validated                          | ✓                                                                                 |                                                                                   |                                                                                   |                                                                                     | ✓                                                                                   | ✓                                                                                   |
| Lot traceable                              | ✓                                                                                 | ✓                                                                                 | ✓                                                                                 | ✓                                                                                   | ✓                                                                                   | ✓                                                                                   |
| <b>Solutions and Packaging</b>             |                                                                                   |                                                                                   |                                                                                   |                                                                                     |                                                                                     |                                                                                     |
| Ready-to-use solution                      | ✓                                                                                 | tablet                                                                            |                                                                                   | ✓                                                                                   | ✓                                                                                   | ✓                                                                                   |
| Concentrate available                      | ✓                                                                                 | tablet                                                                            | ✓                                                                                 |                                                                                     |                                                                                     |                                                                                     |
| Dilution rate for the concentrate solution | 2 oz / gallon                                                                     | 1 tablet / gallon                                                                 | 4 oz / gallon                                                                     |                                                                                     |                                                                                     |                                                                                     |
| Double-bagged bottle/container             | ✓                                                                                 |                                                                                   |                                                                                   | ✓                                                                                   | ✓                                                                                   | ✓                                                                                   |

\*Low levels of VOC materials are in the product: the quat mixture (860 ppm), EDTA (300 ppm) and nonionic surfactant (400 ppm).



**TX<sup>®</sup>650**  
**TexQ<sup>®</sup> Disinfectant**

**GAMMA IRRADIATED**  
Cleaner • Mildewstat  
Fungicide • Sanitizer • Virucide\*  
For Hospitals and Institutional Use

**ACTIVE INGREDIENTS:**  
Octyl Decyl Dimethyl Ammonium Chloride  
Didecyl Dimethyl Ammonium Chloride  
Dioctyl Dimethyl Ammonium Chloride  
Alkyl (C<sub>12</sub>-50%, C<sub>14</sub>-40%, C<sub>16</sub>-10%)  
dimethyl benzyl ammonium chloride  
**INERT INGREDIENTS:**  
**TOTAL**

**KEEP OUT OF REACH OF CHILDREN**  
**CAUTION**  
See Back Panel for Additional Precautions  
and First Aid Statements  
EPA Reg. No. 10324-AB-0000  
Est. No. 48553-AB-0000



# TexQ<sup>®</sup> Disinfectant



## Section Contents

|                                           |   |
|-------------------------------------------|---|
| Why TexQ?.....                            | 4 |
| The TexQ Advantage .....                  | 4 |
| TexQ Products.....                        | 4 |
| Industries & Applications ....            | 4 |
| TexQ Kill Claims .....                    | 5 |
| How TexQ Works.....                       | 6 |
| One-Step Cleaning<br>& Disinfection ..... | 7 |
| Disinfectant Comparison ....              | 8 |
| TexQ Concentrate .....                    | 9 |
| How to Mix .....                          | 9 |

# TexQ Disinfectant

## Why TexQ?

TexQ® Disinfectant is the latest generation of quaternary ammonium compounds (QACs), EPA registered one-step cleaner and disinfectant. It is effective against a broad spectrum of bacteria, viruses and fungi, and inhibits the growth of mold and mildew and their odors when used as directed.

TexQ® Disinfectant is available as: Ready-to-Use spray for small areas (equipment, surfaces) (TX650) and Concentrate solution for large area disinfection (floors, walls) (TX651). Both solutions are 0.2 µm filtered. TexQ® TX650 is gamma-irradiated to a Sterility Assurance Level of 10<sup>-6</sup> according to AAMI Guidelines.

## The TexQ Advantage

- **61 Kill Claims** – Proven efficacy against the most common cleanroom bacteria, viruses and fungi including MRSA, Pseudomonas, H5N1 and HIV-1 at 3 or 10 minutes contact time.
- **Aspergillus Niger Kill Claim** – Most common fungus for clean areas.
- **Complex Formulation** – Excellent cleaning and disinfecting properties for uniform and complete disinfection of surface.
- **One step cleaner/disinfectant** – Cleans and disinfects in one step. Saves additional costs for the cleaner. Saves time and labor. Easy to use.
- **TX650 is Gamma-irradiated** – Compliant with aseptic environment requirements and USP <797>.
- **Free of dyes and fragrance** – No respiratory irritation – safe for the staff, no additional contamination from the vapors.

- **Functional use label on bottle** – Easy documentation and usage control, record the date opened and operator initials.
- **EPA registered** – Kill claims provide assurance of efficacy of disinfection.
- **Hazardous drugs removal** – TX650 is a part of the Hazardous drugs cleaning protocol\* (USP <800>).

## Industries

- Biotechnology
- Hospitals, Pharmacies
- USP <797>, USP <800>
- Medical device manufacturing facilities
- Pharmaceutical manufacturing facilities
- Veterinary clinics and laboratories
- Food processing facilities

## Applications

- Cleaning and disinfecting hard, non-porous surfaces
- Cleaning and disinfecting small surfaces (tables, equipment, isolators, hoods)
- Cleaning and disinfecting large surfaces (floors, walls, ceilings)
- Recommended for use as part of a disinfectant rotation program
- TX650 is recommended for use as part of Texwipe's hazardous drugs cleaning protocol\*

## TexQ Products

| Number |                                                                                     | Description                                                                | Packaging                      |
|--------|-------------------------------------------------------------------------------------|----------------------------------------------------------------------------|--------------------------------|
| TX650  |  | TexQ Disinfectant Ready-to-Use (RTU) Gamma Irradiated 22 oz. Trigger Spray | TX650: 12 bottles/case         |
| TX651  |  | TexQ Disinfectant Concentrate One Gallon                                   | 4 bottles/case beaker included |

\* Hazardous drugs cleaning protocol is available upon request

## Quaternary Ammonium Disinfectant

### TexQ Kill Claims

|                                                                  | TexQ<br>TX650 / TX651<br>Contact Time in<br>Minutes* |
|------------------------------------------------------------------|------------------------------------------------------|
| <b>Bacteria</b>                                                  |                                                      |
| Community Associated Methicillin Resistant Staphylococcus Aureus | 10                                                   |
| Methicillin Resistant Staphylococcus Aureus                      | 10                                                   |
| Burkholderia cepacia                                             | 10                                                   |
| Campylobacter jejuni                                             | 10                                                   |
| Corynebacterium ammoniagenes                                     | 10                                                   |
| Enterobacter aerogenes                                           | 3                                                    |
| Enterobacter cloacae                                             | 10                                                   |
| Enterobacteriaceae w/extended beta lactamase resistance          | 10                                                   |
| Enterococcus faecalis                                            | 10                                                   |
| Enterococcus faecium ( <i>Vancomycin resistant</i> )             | 10                                                   |
| Escherichia coli                                                 | 10                                                   |
| Escherichia coli ( <i>Antibiotic resistant</i> )                 | 10                                                   |
| Escherichia coli O157:H7                                         | 10                                                   |
| Klebsiella pneumoniae                                            | 3                                                    |
| Klebsiella pneumoniae ( <i>Antibiotic resistant</i> )            | 10                                                   |
| Legionella pneumophila                                           | 10                                                   |
| Listeria monocytogenes                                           | 10                                                   |
| Proteus mirabilis                                                | 10                                                   |
| Proteus vulgaris                                                 | 10                                                   |
| Pseudomonas aeruginosa                                           | 10                                                   |
| Pseudomonas aeruginosa ( <i>Antibiotic resistant</i> )           | 10                                                   |
| Salmonella enterica                                              | 10                                                   |
| Salmonella schottmuelleri                                        | 10                                                   |
| Salmonella typhi                                                 | 10                                                   |
| Serratia marcescens                                              | 10                                                   |
| Shigella dysenteriae                                             | 10                                                   |
| Shigella flexneri                                                | 10                                                   |
| Shigella sonnei                                                  | 10                                                   |
| Staphylococcus aureus                                            | 3                                                    |
| Staphylococcus epidermidis ( <i>Antibiotic resistant</i> )       | 10                                                   |
| Streptococcus pyogenes                                           | 10                                                   |
| Vibrio cholerae                                                  | 10                                                   |
| Xanthomonas axonopodis pv. Citri                                 | 10                                                   |
| Xanthomonas campestris pv. Vesicatoria                           | 10                                                   |


|                                                                    | TexQ<br>TX650 / TX651<br>Contact Time in<br>Minutes* |
|--------------------------------------------------------------------|------------------------------------------------------|
| <b>Viruses</b>                                                     |                                                      |
| Avian influenza A Virus ( <i>H5N1</i> )                            | 10                                                   |
| Avian Influenza/Turkey/Wisconsin                                   | 10                                                   |
| Bovine Viral Diarrheal Virus ( <i>BVDV</i> )                       | 10                                                   |
| Canine Coronavirus                                                 | 10                                                   |
| Canine Distemper                                                   | 10                                                   |
| Duck Hepatitis B Virus                                             | 10                                                   |
| Hantavirus                                                         | 10                                                   |
| Hepatitis B virus ( <i>HBV</i> )                                   | 10                                                   |
| Hepatitis C virus ( <i>HCV</i> )                                   | 10                                                   |
| Herpes Simplex Types 1                                             | 10                                                   |
| Herpes Simplex Types 2                                             | 10                                                   |
| HIV-1 ( <i>AIDS virus</i> )                                        | 2                                                    |
| Human Coronavirus                                                  | 10                                                   |
| Infectious Bovine Rhinotracheitis virus ( <i>IBR</i> )             | 10                                                   |
| Influenza Type A / Brazil                                          | 10                                                   |
| Influenza A H1N1 Virus                                             | 10                                                   |
| Newcastle Disease virus                                            | 10                                                   |
| Porcine Respiratory & Reproductive Syndrome Virus ( <i>PRRSV</i> ) | 10                                                   |
| Porcine Rotavirus                                                  | 10                                                   |
| Pseudorabies virus ( <i>Rabies Virus</i> )                         | 10                                                   |
| Respiratory Syncytial ( <i>RSV</i> )                               | 10                                                   |
| Transmissible Gastroenteritis ( <i>TGE</i> )                       | 10                                                   |
| Vaccinia virus ( <i>Pox Virus</i> )                                | 10                                                   |
| <b>Fungi</b>                                                       |                                                      |
| Aspergillus niger                                                  | 10                                                   |
| Candida albicans                                                   | 10                                                   |
| Dactylium dendroides                                               | 10                                                   |
| Trichophyton mentagrophytes ( <i>Athlete's Foot Fungus</i> )       | 10                                                   |



# TexQ Disinfectant

## How Does TexQ Work?

### Disinfection + Cleaning with One Product

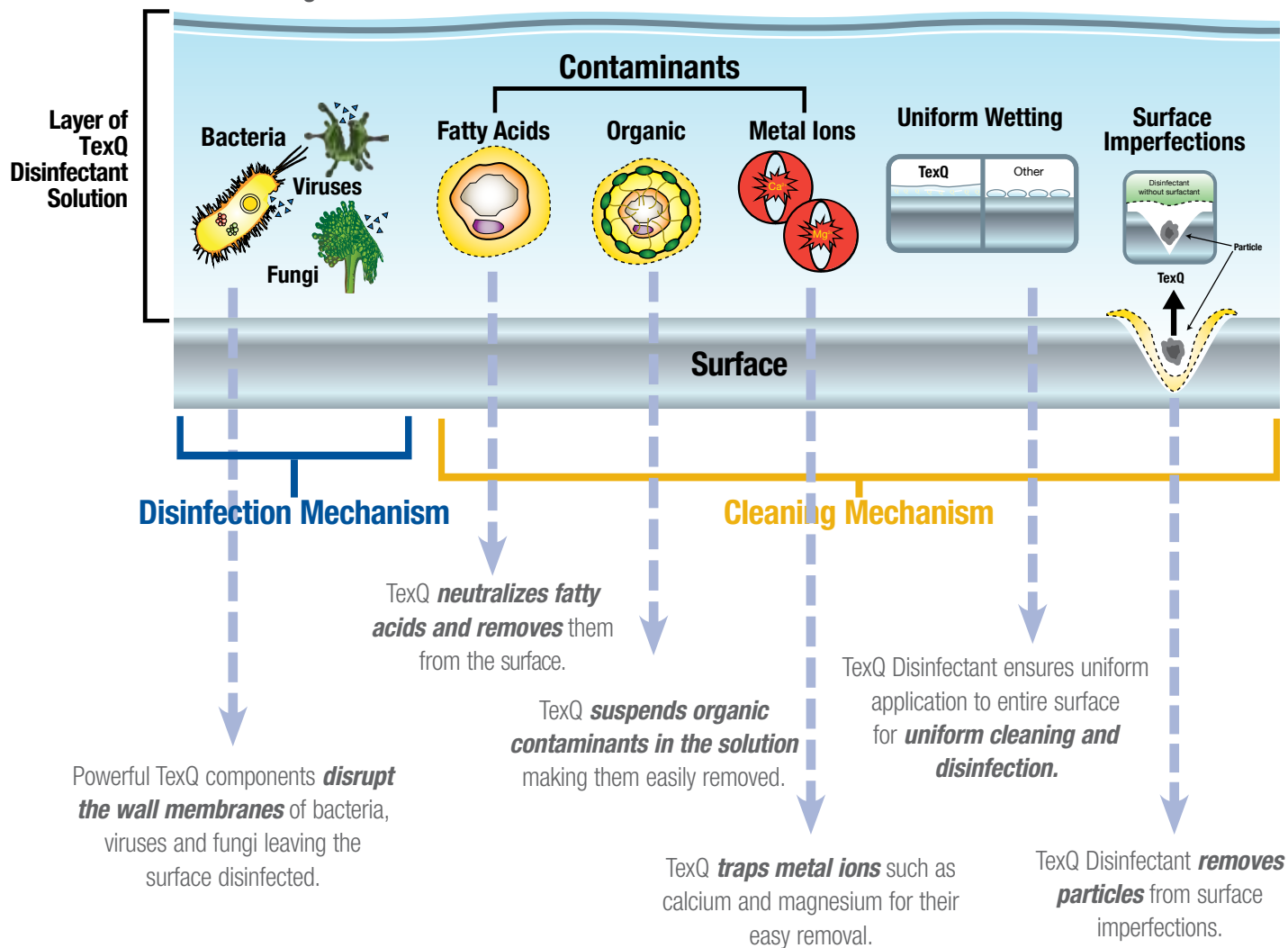

### TexQ Disinfectant is safe on following surfaces:

- Glass, glazed porcelain, glazed ceramic
- Laminated surfaces, Formica®
- Stainless steel, aluminum, metal
- Plexiglass®
- Plastic (such as polycarbonate, polyvinylchloride, polystyrene or polypropylene)
- Vinyl and plastic upholstery
- Sealed granite, sealed marble, sealed limestone, sealed slate, sealed stone
- Sealed terra cotta, sealed terrazzo, vanity tops
- Chrome and vinyl
- Enameled surfaces, painted/finished woodwork
- Washable wallpaper
- External Lenses Vision correction including eyeglasses, protective eyewear, goggles, light lens covers, optical instruments/implements. (Not for use on contact lenses.)

# Quaternary Ammonium Disinfectant

## One-Step Cleaning & Disinfection

### Save a step with TexQ!

The common cleaning and disinfecting practice includes three steps:

1. **Pre-cleaning step** – using a cleaning solution for soil and organic contamination removal.
2. **Disinfectant step** – using a disinfectant solution for antibacterial action.
3. **Residue removal step** – for the disinfectant's left-over removal.

TexQ disinfectant is a one-step disinfectant. That means it **combines the Cleaning and Disinfecting properties** in its complete, chemically balanced formulation. TexQ's **Cleaning Properties** mean that it **removes soil and organic contamination** (by neutralizing fatty acids, suspending organic contaminants and trapping metal ions) from the surface. Its **Disinfecting Properties** provide **antibacterial action** by disrupting the cell walls of bacteria, viruses and fungi.



Just one application of TexQ disinfectant not only disinfects the surface but also removes soil, organic and other contamination combining the “Pre-cleaning step” and the “Disinfection step” in only one easy step\*. The last step remains the “Residue removal” step\*\*. **Having just two steps in the cleaning/disinfection protocol (instead of three) makes the use of TexQ one-step disinfectant solution easier and provides time and labor savings for the operator.**



\* For heavily soiled areas or when visible soil is present, the Pre-cleaning step using a cleaning solution is recommended before any disinfectant application, including one-step disinfectants.

\*\* For non-product contact surfaces (such as walls, floors, ceilings etc) the frequency of the residue removal step may be reduced to once a week/month/two months, according to the approved facility's cleaning procedure.

# TexQ Disinfectant

## TexQ Comparison to Other Disinfectant Classes

|                              | DISINFECTANTS |           |            |                     |
|------------------------------|---------------|-----------|------------|---------------------|
|                              | TexQ          | Bleach    | Phenolics  | Ethanol/Isopropanol |
| Effective pH                 | 8-11          | 10-13     | 1.5 - 12.5 | 6-8                 |
| Cleaning                     | Good          | Poor      | Fair*      | Poor                |
| Need Precleaning Step        | No            | Yes       | No*        | No                  |
| Odor                         | Low           | High      | High       | High                |
| Effectiveness Affected by pH | No            | Yes       | Yes        | No                  |
| Organic Soil Tolerance       | Good          | Poor      | Good*      | Medium              |
| Hard Water Tolerance         | Good          | Good      | Good       | Good                |
| Surface Compatibility        | High          | Medium    | Medium     | High                |
| Corrosiveness                | Low           | High      | Medium     | Low                 |
| Toxicity Category**          | III           | I         | I or II    | IV                  |
| Skin Irritation              | Low           | High      | Medium     | Low                 |
| Respiratory Irritation       | No            | High      | High       | High                |
| Residual Activity            | Yes           | No        | Yes        | No                  |
| Need Rinse                   | Yes           | Yes       | Yes        | No                  |
| Stability / Shelf Life       | Very Good     | Very Poor | Good       | Good                |
| Cost in Use                  | Low           | Low       | Medium     | High                |

\*Depends on product

\*\*EPA Toxicity Categories Require These Warnings:

| Signal Word                           | Category                | Oral Lethal Dose <sup>1</sup>   |
|---------------------------------------|-------------------------|---------------------------------|
| DANGER, POISON (Skull and crossbones) | I Highly toxic          | A few drops to a teaspoonful    |
| WARNING                               | II Moderately toxic     | Over a teaspoonful to one ounce |
| CAUTION                               | III Slightly toxic      | Over one ounce to one pint      |
| CAUTION                               | IV Relatively non-toxic | Over one pint to one pound      |

<sup>1</sup>Based on a 150-pound person

## Shelf Life

For TX650 TexQ Ready-to-use (RTU):

- 2 years (the expiration date is indicated on the product label), even after opening the bottle or spraying from the bottle.

For TX651 TexQ concentrate:

- 2 years (the expiration date is indicated on the product label), even after opening the bottle.
- If the concentrate is diluted to RTU in an open container (or bucket), the solution must be used within 24 hours. If, during use, it gets diluted or visibly dirty, it should be replaced.
- If the concentrate is diluted to RTU and transferred to a spray bottle, it is good for up to 2 years (until the expiration date listed on the original TX651 bottle).

## Disposal

For TX650 and TX651: Dispose of contents and container in accordance with all local, regional, national and international regulations (see Safety Data Sheet, section 13.)

## TexQ Concentrate

**TX651 TexQ Disinfectant** is a concentrated Hospital Use disinfectant effective against a broad-spectrum of bacteria, viruses and fungi, plus eliminates odor-causing bacteria when used as directed. It is available in 1 gallon bottles and features the same properties as the ready-to-use (TX650) spray bottles: 0.2 µm filtered and double-bagged for easy introduction into aseptic areas.

**TexQ Disinfectant Concentrate** formulation meets AOAC Use – Dilution Test Standards for hospital disinfectants and is highly recommended for use in disinfectant rotation programs. At two ounces per gallon of water dilution rate it is proven to be effective against 32 bacteria, 21 viruses and 4 fungi in the presence of organic soil. Treated surfaces must remain wet for 10 minutes. Use this product on large washable, hard, non-porous non-food contact surfaces responsible for cross-contamination such as floors, walls, ceilings, exterior surfaces of the equipment etc.



### How to Mix

Mixing TexQ Disinfectant from TX651 is a simple process: Step 1 – start with 1 gallon of water in the bucket of your choice (we recommend Texwipe Buckets TX7057 Stainless Steel model). Step 2 – Using the included beaker, add 2 ounces of TX651 to the water in the bucket. Step 3 – Mix the 2 ounces of TX651 with the water for 1 minute. TexQ is now mixed and ready to use! *The TexQ concentrate dilution rate is 1:64, or 2 oz per gallon of water.*

#### Step 1



Add 1 gallon of water to bucket.

#### Step 2



Add 2 ounces of TexQ concentrate to the bucket and mix.

#### Step 3



TexQ is ready to use.



**Ideal Buckets for TexQ,** See our [Mops Brochure](#) online (this is a link) at [texwipe.com](http://texwipe.com).





# Bru-Clean TbC™

The Bleach Alternative  
in Tablet Form



## Section Contents

|                                 |    |
|---------------------------------|----|
| Advantages .....                | 12 |
| Applications & Industries ..... | 12 |
| Products .....                  | 12 |
| Comparisons .....               | 13 |
| Kill Claims .....               | 14 |
| How to Mix .....                | 15 |

# Bru-Clean TbC™

Bru-Clean TbC™ is an ideal alternative to liquid bleach that comes in a “just add water” premeasured disinfectant tablet. Bru-Clean TbC™ is designed to provide effective cleaning, deodorizing and disinfection in areas where controlling cross contamination hazards are of the highest concern.

Add 1 Bru-Clean TbC tablet to 1 gallon of water to create a powerful disinfectant with 31 kills claims against a broad array of pathogens including MRSA, E.coli and Salmonella.

The active ingredient, sodium dichloroisocyanurate (NaDCC), is more stable than bleach (sodium hypochlorite) and provides consistent strength produced at point of use, is neutral in pH and is biodegradable.

Bru-Clean TbC meets EPA requirements for hard, inanimate surface disinfectant claims for hospital and medical environments. It is also recommended for use with veterinary, biologics and pharmaceutical manufacturing facilities.

## The Bru-Clean Advantage

- **Convenient premeasured tablets** – Easy to use, transport and store. 270 tablets make 270 gallons of disinfectant.
- **Solution prepared at point-of-use** – Easily prepare a **fresh** solution whenever needed. Consistent concentration of daily prepared solution. Less odor than bleach.
- **Neutral pH** – Much less corrosive to metals than liquid bleach. Less hazardous for users.
- **Shelf life/stability of tablets** – Two years of active ingredient stable concentration in the tablet.
- **EPA registered disinfectant** – Registered kill claims provide the assurance of efficacy of disinfection for 32 microbes.
- **Functional use label on bottle** – Easily document the date opened and operator initials for usage control.

## Applications

- Cleaning and disinfecting hard surfaces, such as stainless steel
- Disinfecting all work areas – equipment, hoods, carts, isolators, floor
- Can easily replace liquid bleach in all applications
- Recommended for use as part of a disinfectant rotation program

## Industries

- Pharmaceutical manufacturing facilities
- Pharmacies
- Hospitals
- Medical device manufacturing facilities
- Biologics and Biotechnology
- Veterinary

## Bru-Clean Products

| Number | Description                         | Packaging                           |
|--------|-------------------------------------|-------------------------------------|
| TX6466 | Bru-Clean TbC™ Disinfectant Tablets | 270 tablets/bottle – 2 bottles/case |



## Liquid Bleach Alternative

### Disinfectant Comparison: Bru-Clean vs Bleach


|                           | Bru-Clean (TX6466)             | Bleach                                        |
|---------------------------|--------------------------------|-----------------------------------------------|
| Effective pH              | 5 - 6                          | 10 - 13                                       |
| Packaging & Delivery      | Bulk packed tablets            | Concentrated solution                         |
| Use convenience           | Made at point-of-use           | Must be stored, diluted, mixed, and filtered  |
| Stability                 | Stable, fresh solution         | Degrades over time (lose 20% in 6 months)*    |
| Shelf Life                | Tablet shelf life of two years | One year                                      |
| Odor                      | Moderate                       | Strong                                        |
| Corrosion potential       | Low                            | High                                          |
| Hazard level: Eye         | Low Irritant                   | High<br>(Severe irritant or may cause damage) |
| Skin                      | Low Irritant                   | High<br>(Severe irritant or may cause damage) |
| Respiratory system        | Irritant                       | High<br>(Severe irritant or may cause damage) |
| DOT Hazard Classification | Not classified as hazardous    | Corrosive, Class 8 at 12% strength            |
| Working concentration     | 937 ppm                        | 2,500 – 5,000 ppm                             |

### Sodium Hypochlorite – Bleach



### Sodium dichloroisocyanurate – NaDCC

Dry tablet, when constituted with water ( $\text{H}_2\text{O}$ )

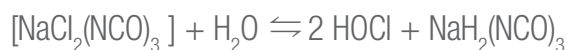

More HOCl is released from NaDCC at pH 5-6 than from sodium hypochlorite (liquid bleach) at pH 9.5

The active ingredient, sodium dichloroisocyanurate (NaDCC), is:

- More stable than bleach (sodium hypochlorite) and provides consistent strength produced at point of use
- Neutral in pH - less corrosive on surfaces than liquid bleach\*
- Biodegradable – safe for the environment

# Bru-Clean TbC™

## Kill Claims

Bru-Clean TbC TX6466  
Contact time (minutes)

| Bacteria                                          |    |
|---------------------------------------------------|----|
| Actinobacillus pleuropneumoniae                   | 10 |
| Brachyspira (Treponema/ Serpuline) hyodysenteriae | 10 |
| Clostridium perfringes USDA                       | 10 |
| Enterococcus faecalis Vancomycin Resistant        | 10 |
| Escherichia coli O157:H7                          | 10 |
| Gumboro disease                                   | 10 |
| Klebsiella pneumoniae                             | 10 |
| Mycobacterium tuberculosis var. bovis             | 10 |
| Pseudomonas aeruginosa                            | 10 |
| Salmonella enterica                               | 10 |
| Staphylococcus aureus                             | 10 |
| Staphylococcus aureus (MRSA)                      | 10 |
| Staphylococcus aureus (GRSA)                      | 10 |
| Staphylococcus epidermidis                        | 10 |
| Streptococcus dysgalactiae                        | 10 |
| Trichophyton mentagrophytes                       | 10 |

Use concentration of 1 tablet per gallon.

Bru-Clean TbC TX6466  
Contact time (minutes)

| Viruses                                  |    |
|------------------------------------------|----|
| Adenovirus Type 2                        | 10 |
| Avian Influenza Type A                   | 10 |
| Canine distemper virus                   | 10 |
| Canine Parvovirus                        | 10 |
| Feline calicivirus                       | 10 |
| Hepatitis A virus (HAV)                  | 10 |
| Herpes Simplex Type 1                    | 10 |
| HIV-1 (AIDS virus)                       | 10 |
| Infectious canine hepatitis              | 10 |
| Newcastle disease virus                  | 10 |
| Norovirus                                | 10 |
| Poliovirus type 1                        | 10 |
| Porcine parvovirus                       | 10 |
| Pseudorabies                             | 10 |
| Runting & Stunging virus (tenosynovitis) | 10 |
| Fungi                                    |    |
| Trichophyton mentagrophytes              | 10 |

## Liquid Bleach Alternative

### How to Mix

According to the AOAC Use Dilution Test method, this product is bactericidal, virucidal and fungicidal with 32 registered kill claims. Treated surfaces must remain wet for 10 minutes. Concentration of available chlorine – 937 ppm (1 tablet per gallon of water dilution rate).



**Step 1** Add 1 gallon of water to bucket.



**Step 2** Add one tablet then wait 2 to 3 minutes while it fizzes & dissolves.



**Step 3** Mix and your disinfectant solution is ready to use!



**Ideal Buckets for Bru-Clean TbC,**

See our [Mops Brochure](#) online (this is a link) at [texwipe.com](http://texwipe.com).



# TexCide™



## Section Contents

Why Texcide? .....18

Features & Benefits .....18

Industries & Applications ...18

TexCide Products .....18

TexCide Kill Claims .....19

How does TexCide  
kill spores? .....20

Texwipe's Cleaning Rotation  
Recommendations .....21

# TexCide™ TX690

## Why TexCide?

TexCide™ TX690 is an effective EPA registered 5.9% peroxyacetic acid/27.3% hydrogen peroxide-based sporicidal disinfectant with fast efficacy against *Clostridium difficile* (C. diff) spores.

TexCide is a concentrated, broad-spectrum, one-step disinfectant cleaner and deodorizer effective against spores, bacteria, viruses and fungi just within a two (2) minute contact time. The 8 oz. (250 ml) measuring beaker makes the dilution process easier.

- **Fragrance free.**
- **Non-flammable.**
- **Measuring beaker included** – makes the dilution process easier.
- **Lot traceable.**

*\*at the dilution rate 4 oz. per gallon of water (see the kill claims on reverse)*

*\*\*if no visible soil is present*

## Features & Benefits

- **Fast sporicidal action** – Kills *Clostridium difficile* spores in 2 minutes.\*
- **Bactericidal, virucidal, and fungicidal** with a contact time of 2-5 minutes.\*
- **One-step disinfectant and cleaner** – cleans, disinfects and deodorizes in one labor saving step.
- **No pre-cleaning step is required \*\*** – effective in the presence of 5% organic serum load. Saves additional costs incurred from a pre-cleaning step (cleaning ingredient, time, etc.).
- **EPA registered.**
- **Compatible with stainless steel** and other 11 surfaces (see next page).
- **No special water requirements for dilution** – effective in water up to 400 ppm hardness.
- **No rinsing required** when used on floors, walls and ceilings.
- **Convenient to use** – 16 oz. bottle of concentrate makes 4 gallons of use solution.

## Industries

- Pharmaceutical
- Biotechnology
- Hospitals and Healthcare facilities
- Medical Device
- Laboratories
- Pharmacies and Compounding Pharmacies
- Veterinary clinics

## Applications

- Cleaning and disinfecting hard, non-porous surfaces
- Cleaning and disinfecting small surfaces (tables, equipment, isolators, hoods)
- Cleaning and disinfecting large surfaces (floors, walls, ceilings)
- Recommended for use as part of a disinfectant rotation program as the sporicidal agent

## TexQ Products

| Number | Description          | Packaging                                                     |
|--------|----------------------|---------------------------------------------------------------|
| TX690  | TexCide™ Concentrate | 16 oz (473 ml) concentrate, 12 bottles/case (beaker included) |



## TexCide Kill Claims

|                                                                          | TexCide TX690<br>Contact Time<br>in Minutes* |                                                                           | TexCide TX690<br>Contact Time<br>in Minutes* |
|--------------------------------------------------------------------------|----------------------------------------------|---------------------------------------------------------------------------|----------------------------------------------|
| <b>Sporicidal kill claims</b>                                            |                                              | <b>Virucidal kill claims</b>                                              |                                              |
| Clostridium difficile                                                    | 2                                            | Adenovirus Type 5                                                         | 2                                            |
| <b>Bactericidal kill claims</b>                                          |                                              | Canine Parvovirus (CPV)                                                   | 5                                            |
| Acinetobacter baumannii                                                  | 2                                            | Hepatitis B Virus (HBV)                                                   | 5                                            |
| Bordetella pertussis                                                     | 2                                            | Hepatitis C Virus (HCV)                                                   | 5                                            |
| Enterococcus faecalis Vancomycin Resistant (VRE)                         | 2                                            | Herpes Simplex Virus Type 1                                               | 2                                            |
| Escherichia coli                                                         | 2                                            | Herpes Simplex Virus Type 2                                               | 2                                            |
| Escherichia coli with beta-lactamase resistance (ESBL)                   | 2                                            | Human Immunodeficiency Virus Type 1 (HIV-1)                               | 2                                            |
| Klebsiella pneumoniae                                                    | 2                                            | Influenza A Virus                                                         | 2                                            |
| Klebsiella pneumoniae Carbapenem Resistant                               | 2                                            | Murine Norovirus (MNV-1)                                                  | 2                                            |
| Proteus mirabilis                                                        | 2                                            | Norovirus                                                                 | 2                                            |
| Pseudomonas aeruginosa                                                   | 2                                            | Respiratory Syncytial Virus (RSV)                                         | 2                                            |
| Salmonella enterica                                                      | 2                                            | Rhinovirus Type 37                                                        | 2                                            |
| Staphylococcus aureus                                                    | 2                                            | Rotavirus                                                                 | 2                                            |
| Staphylococcus aureus Methicillin Resistant (MRSA)                       | 2                                            | Vaccinia Virus                                                            | 2                                            |
| Staphylococcus aureus Community Acquired Methicillin Resistant (CA-MRSA) | 2                                            | <b>Fungicidal kill claims</b>                                             |                                              |
| Staphylococcus aureus Vancomycin Intermediate Resistant (VISA)           | 2                                            | Candida albicans                                                          | 2                                            |
| Streptococcus pneumoniae                                                 | 2                                            | Trichophyton mentagrophytes (Athlete's foot fungus) (a cause of Ringworm) | 2                                            |
| Streptococcus pyogenes                                                   | 2                                            | TOTAL                                                                     | 33                                           |

\*Tested according to the AOAC Use Dilution method on hard, non-porous surfaces, at 4 ounces per gallon of 400 ppm hard water (1844 ppm active PAA) in the presence of 5% organic serum.

## Typical Properties

- Water solubility Complete
- Physical form Liquid
- Color Colorless
- pH (25°C) 2.65
- Dilution recommendation — 4 oz. per gallon of water

## Surface compatibility


- Finished floors
- Enameled surfaces
- Desks
- Glass
- Glazed porcelain
- Medical equipment surfaces
- Plastic & painted surfaces
- Stainless steel
- Vinyl
- Aluminum
- Chrome
- Examination tables
- Glazed ceramic
- Laminated surfaces
- Workstations

## Not recommended for use on

- Copper
- Brass
- Granite
- Marble
- Zinc
- Unsealed/uncoated marble or terrazzo floors

# TexCide™ TX690

## How does TexCide kill spores?

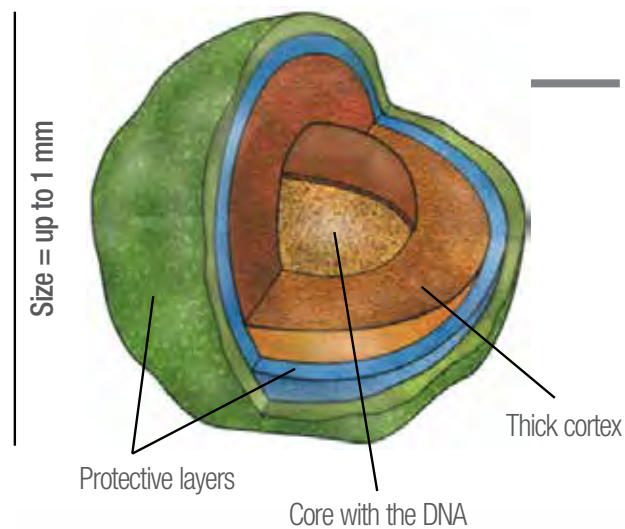

**TexCide kills a spore in just 2 minutes!**

### Sporicidal action:

- TexCide's active ingredients kill the spore by disrupting its protective layers, thick cortex and the spore core.

### Bactericidal, virucidal, fungicidal action:

- TexCide kills other bacteria, viruses and fungi in just 2 minutes!



UV light



Heat >105°F  
(>40°C)



Common chemicals

A thick cortex and protective layers protect the spore from UV light, heat and chemicals

| PRODUCT                                    | TexCide TX690                                         | Spor-Klenz Concentrate Sterilant                    | Spor-Klenz RTU                                      | Decon Spore 200 Plus                                  | SporGon                                                | Peridox RTU                                           |
|--------------------------------------------|-------------------------------------------------------|-----------------------------------------------------|-----------------------------------------------------|-------------------------------------------------------|--------------------------------------------------------|-------------------------------------------------------|
| RTU or Concentrate                         | Concentrate                                           | Concentrate                                         | RTU                                                 | Concentrate                                           | RTU                                                    | RTU                                                   |
| <b>Sporicidal contact time, in minutes</b> | <b>2</b>                                              | <b>No spore kill claim</b>                          | <b>30</b>                                           | <b>3</b>                                              | <b>180</b>                                             | <b>3</b>                                              |
| Contact time, min.                         |                                                       |                                                     |                                                     |                                                       |                                                        |                                                       |
| Bactericidal                               | 2                                                     | 10                                                  | 10                                                  | 10                                                    | 3-15                                                   | 2-5                                                   |
| Virucidal                                  | 2                                                     | No kill claims                                      | 10                                                  | 10                                                    | 5                                                      | 2-3                                                   |
| Fungicidal                                 | 2                                                     | No kill claims                                      | 5                                                   | No kill claims                                        | 5                                                      | 1                                                     |
| <b>Total number of kill claims</b>         | <b>33</b>                                             | <b>2</b>                                            | <b>15</b>                                           | <b>10</b>                                             | <b>10</b>                                              | <b>51</b>                                             |
| Product formulation                        | Hydrogen Peroxide - 27.3%<br>Peroxyacetic acid - 5.9% | Hydrogen Peroxide - 22%<br>Peroxyacetic acid - 4.5% | Hydrogen Peroxide - 1%<br>Peroxyacetic acid - 0.08% | Hydrogen Peroxide - 27.5%<br>Peroxyacetic acid - 5.8% | Hydrogen Peroxide - 7.35%<br>Peroxyacetic acid - 0.23% | Hydrogen Peroxide - 4.4%<br>Peroxyacetic acid - 0.23% |
| Available sizes                            | 16 oz (12/case)                                       | 32 oz (4/case)                                      | 29 oz (4/case)                                      | various                                               | 1 gallon (4/case)                                      | 32 oz (6/case)                                        |



### TexCide:

- has the **shortest sporicidal contact time** and one of the largest total number of kill claims vs the compared products
- the bactericidal, virucidal and fungicidal kill times are the same as the sporicidal kill time (2 min) — for **more efficient use of disinfection time**
- concentrated solution is more **cost effective** than the RTU

### Products having no EPA-registered sporicidal kill claims:

- Spor-Klenz Concentrate (Steris)
- Accel Tb, Accel Concentrate (Contec)
- Steri-Perox (Veltek)
- Clorox Healthcare® Hydrogen Peroxide
- Cleaner Disinfectant (Clorox)
- Hypo-Chlor (Veltek)
- Sporicidin (Contec)

# Texwipe's Cleaning Rotation Recommendations
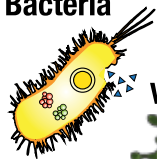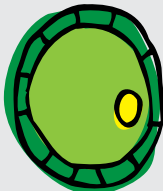
| Microbe to be killed                                                                                                                                                                                                                                                                                                                                           | Disinfectants to be used                                                                                                                                                                                                                                                                                                 | How often to use                                                                                                                                                              | Remove disinfectant residues                                                                                                                                                                                                                                             |
|----------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|--------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|-------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|--------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
| <b>Bacteria</b><br><br><b>Viruses</b><br><br><b>Fungi</b><br><br><b>Vegetative Bacteria, viruses, fungi</b> | <b>Bactericidal Disinfectants</b> <ul style="list-style-type: none"> <li>Quaternary Ammonium Compounds (QAC)<br/><i>TexQ disinfectant</i></li> <li>and</li> <li>Chlorine compound<br/>&lt;5000 ppm<br/><i>Bru-Clean TbC</i></li> </ul>  | <b>Use daily</b> (for example, at the beginning and/or at the end of the shift);<br><br><b>Change weekly</b> (for example, 1st week – use TexQ, 2nd week – use Bru-Clean TbC) | <br><b>TexP</b><br>Hydrogen Peroxide Solution<br><br><br><b>70% IPA</b><br>sterile or non-sterile |
| <br><b>Bacterial Spores (C.Diff)</b>                                                                                                                                                                                                                                         | <b>Sporicidal Disinfectants</b> <ul style="list-style-type: none"> <li>Hydrogen peroxide/peracetic acid<br/><i>TexCide</i></li> </ul>                                                                                                  | Use once a week, or every 2 weeks (for example, on Fridays, at the end of the last shift)                                                                                     | <br><b>70% IPA</b><br>sterile or non-sterile                                                                                                                                         |


## Dilution recommendations:

- 4 oz. of concentrate per gallon of water.

Use a full 16 oz bottle to make a 4 gallon bucket of use solution (without using the beaker):

1. Add 4 gallons of water to bucket
2. Empty a 16 oz bottle in the bucket and mix
3. Your sporicidal solution is ready



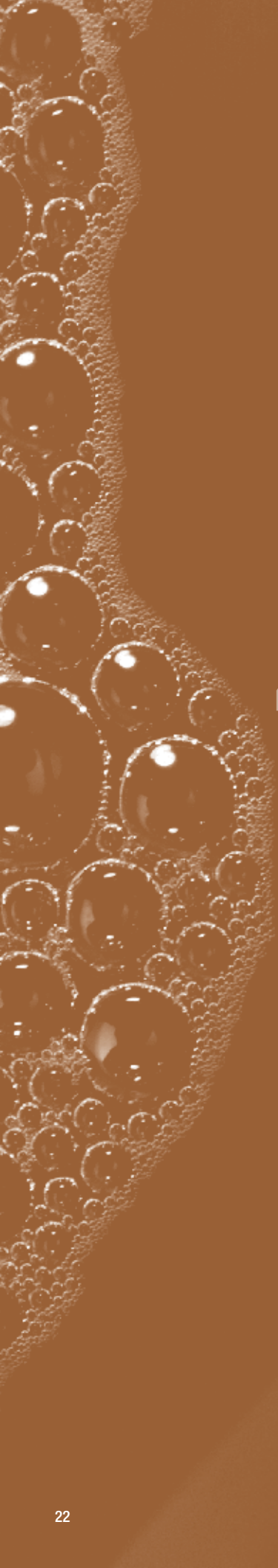

4%

7.5%

# TexP™

## Non-Sterile Hydrogen Peroxide



### Section Contents

|                           |    |
|---------------------------|----|
| TexP .....                | 24 |
| Features & Benefits ..... | 24 |
| TexP Products .....       | 24 |
| Industries .....          | 25 |
| Applications .....        | 25 |

# TexP

## Hydrogen Peroxide, Non-Sterile

Texwipe's® TexP (hydrogen peroxide) cleaning solutions are developed for use with cleaning rotation cycles to maintain a clean environment. TexP offers no VOCs and is nonflammable, making it safer for your personnel and regulatory compliance.

TexP leaves no residue, making it ideal for residue removal with disinfectants such as TexQ, Bru-Clean TbC and other formulations. It offers a low toxicity profile and is noncorrosive on most surfaces. TexP is completely biodegradable, it completely decomposes into water and oxygen.

TexP is offered in two ready-to-use concentrations – 4% and 7.5% – and both are available in 16 oz. spray (suitable for small areas) and 1 gallon (ideal for large volume applications) bottles. All TexP products are individually lot-coded for ease of traceability and quality control.

Labeled in three languages (English, Spanish, French).

## TexP Advantages

- **Convenient ready-to-use solutions** require no mixing or activating.
- **Nonflammable** – may be used in high flammability risk environments.
- **No VOCs** – reduces VOC regulatory and compliance concerns.
- **No residue**, may be used for disinfectant removal.
- **Safe for use** – low toxicity profile for staff and noncorrosive on most common surfaces.
- **Biodegradable** – completely decomposing into water and oxygen.

## TexP Products

| Solution               | Number |                                                                                     | Description                                                    | Packaging           |
|------------------------|--------|-------------------------------------------------------------------------------------|----------------------------------------------------------------|---------------------|
| 4% hydrogen peroxide   | TX684  |  | 4% hydrogen peroxide<br>Ready-to-use<br>Non-sterile solution   | 16 oz. spray bottle |
|                        | TX684G |  | 4% hydrogen peroxide<br>Ready-to-use<br>Non-sterile solution   | 1 gallon bottle     |
| 7.5% hydrogen peroxide | TX687  |  | 7.5% hydrogen peroxide<br>Ready-to-use<br>Non-sterile solution | 16 oz. spray bottle |
|                        | TX687G |  | 7.5% hydrogen peroxide<br>Ready-to-use<br>Non-sterile solution | 1 gallon bottle     |

## 4% and 7.5% Hydrogen Peroxide

### Industries

- Pharmaceuticals
- Biologics and Biotechnology
- Medical Device
- Compounding Pharmacy
- Cosmetics
- Health Care
- Veterinary
- Food and Beverage
- Nutraceutical
- Microelectronics
- Semiconductor

### Applications

- Cleaning small (tables, equipment) surfaces
- Cleaning large (floors, walls, ceilings) surfaces
- Removing residues after disinfectant application





# IPA

## Sterile & Non-Sterile Isopropyl Alcohol



### Section Contents

|                    |    |
|--------------------|----|
| 70% IPA .....      | 28 |
| 100% IPA .....     | 28 |
| IPA Products ..... | 29 |

# IPA Solutions

---

## 70% and 100%

Available in both Sterile and Non-Sterile form, Texwipe's Isopropyl Alcohol Solutions are 0.2 micron filtered and filled into cleaned, double-bagged containers. Each bottle is lot coded with an expiration date for easy record keeping.

The trigger spray bottles are fully assembled and ready to use - no unpacking, handling or attaching a separate sprayhead is required. Choose a stream delivery for precise application or coarse spray for larger area application. Gallon containers are also available.

## Features & Benefits

- 0.2 µm filtered Isopropanol (Isopropyl Alcohol) – removes many surface contaminants. (70% is USP-grade and 100% is semiconductor-grade)
- Evaporates and leaves extremely low residue
- Adjustable trigger spray – allows the liquid to be dispensed as either a jet or spray.
- Double-bagged for easy introduction into clean environment
- Meets FDA guideline for medical device bacterial endotoxin limit of not more than 20 EU/device
- Fully lot traceable
- Eliminates mixing, filtration, processing and QC expense of in-house production of sterile IPA

### Sterility Assurance

- Gamma-irradiated to a sterility level of 10<sup>-6</sup> with independent QC audits to ensure sterility
- Each lot tested for endotoxins
- Lot traceable
- Lot-specific information on each shipment simplifies record keeping:
  - » Certificate of Processing confirming radiation dosage
  - » Certificates of Compliance & Analysis

## Applications

- Surface cleaning
- Residue removal after disinfectant application
- Use on gloved hands
- Wipe down for pass through to controlled environments
- Ideal for use with cleanroom wipers

---

## 70% IPA Solutions

Texwipe's 70% Isopropanol contains 70% by volume USP-grade Isopropanol (Isopropyl Alcohol) and 30% USP-purified water.

### Industries

- Pharmaceuticals
- Biologics
- Medical Device
- Compounding Pharmacy

---

## 100% IPA Solutions

Non-Sterile 100% Isopropyl Alcohol is made with semiconductor-grade Isopropanol.

### Industries

- Microelectronics

## IPA Products

| IPA %                         | Sterile | Number           |                                                                                     | Description                                                                        | Packaging                                                 |
|-------------------------------|---------|------------------|-------------------------------------------------------------------------------------|------------------------------------------------------------------------------------|-----------------------------------------------------------|
| <b>70% Isopropyl Alcohol</b>  |         |                  |                                                                                     |                                                                                    |                                                           |
| 70%                           | ●       | TX8270           |    | 70% Isopropyl Alcohol<br>Trigger Spray<br>8 fluid ounces (237 mL)<br>Sterile       | 12 polybottles per case                                   |
| 70%                           | ●       | TX3270<br>TX3274 |    | 70% Isopropyl Alcohol<br>Trigger Spray<br>16 fluid ounces (473 mL)<br>Sterile      | TX3270: 12 bottles per case<br>TX3274: 4 bottles per case |
| 70%                           |         | TX167            |    | 70% Isopropyl Alcohol<br>Trigger Spray<br>16 fluid ounces (473 mL)<br>Non-Sterile  | 12 bottles per case                                       |
| 70%                           | ●       | TX3273           |   | 70% Isopropyl Alcohol<br>Trigger Spray<br>32 oz. spray bottle (946 mL)<br>Sterile  | 12 bottles per case                                       |
| 70%                           | ●       | TX3290           |  | 70% Isopropyl Alcohol<br>1 gallon (3.8 liters)<br>Sterile                          | 4 bottles per case                                        |
| 70%                           |         | TX117            |  | 70% Isopropyl Alcohol<br>1 gallon (3.8 liters)<br>Non-Sterile                      | 4 bottles per case                                        |
| <b>100% Isopropyl Alcohol</b> |         |                  |                                                                                     |                                                                                    |                                                           |
| 100%                          |         | TX161            |  | 100% Isopropyl Alcohol<br>1 gallon (3.8 liters)<br>Non-Sterile                     | 4 bottles per case                                        |
| 100%                          |         | TX111            |  | 100% Isopropyl Alcohol<br>Trigger Spray<br>16 fluid ounces (473 mL)<br>Non-Sterile | 12 bottles per case                                       |

# ETOH

## Denatured 70% Ethanol



# 70% Denatured Ethanol

Sterile, ready-to-use 70% denatured ethanol (EtOH) solution designed especially for cleanroom use. Filtered to 0.2µm, individually double-bagged and gamma-irradiated to ensure sterility. Made using USP components and packaged in a ISO Class 5 cleanroom.

## Features & Benefits

- **Filled in an ISO Class 5 environment;** 0.2µm filtered
- **Individually double bagged and gamma irradiated** to a Sterility Assurance Level of 10<sup>-6</sup> with independent QC audits to ensure sterility
- **Evaporates** leaving extremely low residue
- **Adjustable trigger spray** which allows the liquid to be dispensed as either a jet or spray
- **Fully lot traceable.** Each lot tested for endotoxins
- **Each shipment is supported by documentation:** Certificate of Irradiation, Certificate of Compliance and Analysis
- **MSDS and Sterility Validation** documents available

## Industries

- Microelectronics
- Semiconductors
- Biologics
- Medical Device
- Pharmaceuticals
- Compounding Pharmacies

## Applications

- Surface cleaning and residue removal purposes
- Use on gloved hands in sterile suites
- Wipe down for pass through to controlled environments
- Ideally suited for use with sterile cleanroom wipers

## Sterility Assurance

- Gamma-irradiated to a sterility level of 10<sup>-6</sup> with independent QC audits to ensure sterility
- Each lot tested for endotoxins
- Lot traceable
- Lot-specific information on each shipment simplifies record keeping:
  - » Certificate of Processing confirming radiation dosage
  - » Certificates of Compliance & Analysis

## 70% Ethanol Products

| Number         | Description                                                                                                       | Packaging                                                      |
|----------------|-------------------------------------------------------------------------------------------------------------------|----------------------------------------------------------------|
| <b>Sterile</b> |                                                                                                                   |                                                                |
| TX3267         |  Sterile 70% Denatured Ethanol | 16 oz. trigger spray bottle (473 mL) / 12 polybottles per case |
| TX3265         |  Sterile 70% Denatured Ethanol | 32 oz. trigger spray bottle (946 mL) / 12 polybottles per case |



# CrushTube<sup>TM</sup> Swab

91% IPA / 9%DIW

# CrushTube™ Swab

Texwipe's CrushTube system includes a 100% polyester (hydroentangled) nonwoven material head attached to an internal vial containing 91% IPA / 9% DIW solution, which is enclosed in a protective casing.

When the internal vial is crushed, the head becomes saturated with the 91% IPA / 9% DIW solution for point-of-use application.

Individually packaged in cleanroom compatible material. Lot coded for traceability and quality control.

## Features & Benefits

- **IPA solution is separated from head until activated**
- **Gently crush** the vial to release the IPA and saturate the brush/tip
- **Precision spot cleaning**, no extra container of solvent needed
- **Individually packaged to minimize storage need**, easy to transport and use

## Applications

- Solvent cleaning and maintaining of ion emitter tips
- Cleaning of grooves, tracks, slots and other small spaces
- Removing adhesive buildup
- Solvent cleaning sensitive surfaces such as optical assemblies

## Industries

- Automotive
- Biologics
- Food Manufacturing
- Microelectronics
- Pharmaceuticals
- Semiconductor

## CrushTube Products

| Number |                                                                                     | Description | Packaging                   |
|--------|-------------------------------------------------------------------------------------|-------------|-----------------------------|
| TX726  |  | CrushTube™  | 50 swabs/box; 10 boxes/case |

**For More Swab Solutions,** See our [Swabs Brochure](#) online (this is a link) at [texwipe.com](http://texwipe.com).



# APPENDIX

## Section Contents

Disinfectant  
Classification . . . . . 35

Disinfectant + Wiper  
Compatibility . . . . . 36

Pre-Wetted Wipers. . . . . 37

ISO 5 Sample  
Cleaning Frequency. . . . . 38



# Disinfectants Classification

|                                                   | Isopropyl Alcohol                                                                                                                                                                                                                                                                                                                                                        | Chlorine Compounds, Bleach Sodium Hypochlorite 5.25% (bleach concentrate)                                                                                                                                                                                                                                                                                                                                                                                                       | Phenolics                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                      | Quaternary Ammonium Compounds (QACs)                                                                                                                                                                                                                                                                                                                                                                                                                    | Oxidizing Disinfectants, Hydrogen peroxide                                                                                                                                                                                                                                                                                                                                                                                           |
|---------------------------------------------------|--------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|---------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|--------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|---------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|--------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
| <b>Description</b>                                | Variable activity against some bacterial and fungal species. Bactericidal disinfectant.<br>70% IPA is proved to be the most effective concentration.                                                                                                                                                                                                                     | Bactericidal (kills bacteria, viruses and fungi) at <5000ppm Sodium Hypochlorite.<br>Sporicidal (kills spores) at >5000ppm Sodium Hypochlorite.                                                                                                                                                                                                                                                                                                                                 | Bactericidal disinfectant (kills bacteria, viruses, fungi), Tuberculocidal.                                                                                                                                                                                                                                                                                                                                                                                                                                                    | Bactericidal disinfectant (kills bacteria, viruses and fungi). Some products are Tuberculocidal.                                                                                                                                                                                                                                                                                                                                                        | This group includes oxygen-releasing compounds (peroxygens) such as peracetic acid and hydrogen peroxide.<br>Bactericidal (kills bacteria, viruses, fungi), Tuberculocidal. Sporicidal (kills spores).                                                                                                                                                                                                                               |
| <b>Pre-cleaning Needed</b>                        | Surfaces must be pre-cleaned.                                                                                                                                                                                                                                                                                                                                            | Surfaces must be pre-cleaned.                                                                                                                                                                                                                                                                                                                                                                                                                                                   | Surfaces must be pre-cleaned. Some products are registered as one step disinfectant cleaners.                                                                                                                                                                                                                                                                                                                                                                                                                                  | Product specific. Some products registered as one-step disinfectant cleaners.                                                                                                                                                                                                                                                                                                                                                                           | Product specific. Surfaces must be pre-cleaned, depending on formulation.                                                                                                                                                                                                                                                                                                                                                            |
| <b>Advantages</b>                                 | <ul style="list-style-type: none"> <li>• Quick evaporation</li> <li>• Removes many surface contaminants</li> <li>• Removes residual disinfectant</li> <li>• Leaves extremely low residue</li> <li>• Can be a good general use disinfectant</li> <li>• Compatibility combined with other disinfectants (quaternaries, phenolics)</li> <li>• No rinse required</li> </ul>  | <ul style="list-style-type: none"> <li>• The same product can be used for routine and special event tasks, by changing the concentration</li> <li>• Relatively quick microbial kill</li> <li>• May be used on food preparation surfaces requiring a surface rinse depending on bleach concentration</li> <li>• Can be Tuberculocidal and Sporicidal with increased concentration</li> <li>• Some products are Tuberculocidal</li> </ul>                                         | <ul style="list-style-type: none"> <li>• Mostly presented in concentrate formulations, need to be mixed to make the ready-to-use solution</li> <li>• Some products are Tuberculocidal</li> <li>• Effective over large pH range</li> <li>• Some products are one-step disinfectants cleaners</li> </ul>                                                                                                                                                                                                                         | <ul style="list-style-type: none"> <li>• One-step formulations contain a detergent to help loosen soil, no pre-clean step needed</li> <li>• Colorless, odorless (but act as deodorizers)</li> <li>• Less corrosive</li> <li>• May be used on food preparation surfaces (need rinse)</li> <li>• Effective at temperatures up to 212°F</li> </ul>                                                                                                         | <ul style="list-style-type: none"> <li>• Hydrogen Peroxide is non-corrosive in diluted form but is corrosive in combination with peracetic acid</li> <li>• No rinsing required</li> <li>• Some products are odorless</li> <li>• Clear and colorless, thereby avoiding surface staining</li> <li>• Fast, broad spectrum activity, Sporicidal</li> <li>• Can be safer for personnel (less toxic) depending on concentration</li> </ul> |
| <b>Disadvantages</b>                              | <ul style="list-style-type: none"> <li>• Poor cleaner (does not contain detergents)</li> <li>• Limited contact time, not sufficient for broad range killing, evaporates quickly</li> <li>• VOC emissions</li> <li>• Flammable, not to be used near a flame</li> <li>• Not active against certain types of viruses</li> <li>• Low toxicity but an eye irritant</li> </ul> | <ul style="list-style-type: none"> <li>• Toxic. May damage floor finishes, carpets, clothing and other fibers when used in higher concentrations</li> <li>• Has an unpleasant odor</li> <li>• Must be stored separately from ammonia and flammable products</li> <li>• Rinsing is required</li> <li>• Corrodes metals such as stainless, aluminum</li> <li>• Increase in alkalinity decreases bactericidal properties</li> <li>• Eye, skin, and respiratory irritant</li> </ul> | <ul style="list-style-type: none"> <li>• Considered a persistent bio accumulative toxin by EPA</li> <li>• Disposal restrictions in some states. Check state and local regulations</li> <li>• Not for use on food or food utensils</li> <li>• May damage floor finishes and other surfaces</li> <li>• Unpleasant odor</li> <li>• Effectiveness reduced by alkaline pH or natural soap</li> <li>• Prolonged contact deteriorates rubber</li> <li>• Can cause skin and eye irritation</li> <li>• Corrosive &amp; toxic</li> </ul> | <ul style="list-style-type: none"> <li>• Ineffective against bacterial spores, TBC, some viruses</li> <li>• Effectiveness influenced by hard water</li> <li>• RTU formulations are non-irritating to skin but avoid skin or eye contact; toxic</li> <li>• Neutralized by anionic soap (common) and effectiveness reduced by organic material</li> <li>• Pre-rinse may be required when rotating disinfectants</li> <li>• Rinsing is required</li> </ul> | <ul style="list-style-type: none"> <li>• Rinsing is required where direct skin or oral contact can occur</li> <li>• Corrosive to soft metals</li> <li>• Pre-cleaning step is required</li> <li>• Temperature and light sensitive</li> <li>• Pungent odor (vinegar)</li> <li>• Pure Hydrogen Peroxide formulations do not require rinse</li> </ul>                                                                                    |
| <b>CDC Disinfection Level</b>                     | Intermediate                                                                                                                                                                                                                                                                                                                                                             | Intermediate Level Disinfectant                                                                                                                                                                                                                                                                                                                                                                                                                                                 | Some are intermediate some are low level noted on label                                                                                                                                                                                                                                                                                                                                                                                                                                                                        | Low Level Disinfectant                                                                                                                                                                                                                                                                                                                                                                                                                                  | Product Specific. Low, Intermediate or High Level Disinfectant (depends on concentration and exposure time).                                                                                                                                                                                                                                                                                                                         |
| <b>EPA Toxicity Category</b><br>(See chart below) | Category IV                                                                                                                                                                                                                                                                                                                                                              | Category I                                                                                                                                                                                                                                                                                                                                                                                                                                                                      | Category I or II                                                                                                                                                                                                                                                                                                                                                                                                                                                                                                               | Category III                                                                                                                                                                                                                                                                                                                                                                                                                                            | Category III or IV, product specific.                                                                                                                                                                                                                                                                                                                                                                                                |
| <b>Storage</b>                                    | Stable in storage. Keep away from oxidizing agents, heat and flames.                                                                                                                                                                                                                                                                                                     | If used for disinfecting purposes, bleach should not be stored longer than 3 months. When mixed with water the solution is only effective as a disinfectant for 24 hours. The available chlorine level (NaOCl) must be monitored.                                                                                                                                                                                                                                               | Stable in storage. Flammable if in aerosol form.                                                                                                                                                                                                                                                                                                                                                                                                                                                                               | Stable in storage.                                                                                                                                                                                                                                                                                                                                                                                                                                      | Stable in storage. Two year shelf life is available depending on concentration and formulation. Keep away from heat and light.                                                                                                                                                                                                                                                                                                       |

\*\*EPA Toxicity Categories Require These Warnings:

| Signal Word                           | Category                | Oral Lethal Dose <sup>1</sup>   |
|---------------------------------------|-------------------------|---------------------------------|
| DANGER, POISON (Skull and crossbones) | I Highly toxic          | A few drops to a teaspoonful    |
| WARNING                               | II Moderately toxic     | Over a teaspoonful to one ounce |
| CAUTION                               | III Slightly toxic      | Over one ounce to one pint      |
| CAUTION                               | IV Relatively non-toxic | Over one pint to one pound      |

<sup>1</sup>Based on a 150-pound person

# Disinfectant + Wiper Product Compatibility

| Material/Fabric             | Texwipe Wiper Products                                                                                                                                                                                                                                     | Texwipe Mop Products                                                                                      | Disinfectants |               |         | Cleaners |     |         |
|-----------------------------|------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|-----------------------------------------------------------------------------------------------------------|---------------|---------------|---------|----------|-----|---------|
|                             |                                                                                                                                                                                                                                                            |                                                                                                           | TexQ          | Bru-Clean TbC | Texcide | TexP     | IPA | Ethanol |
| Polyester/<br>Cellulose     | TX622, 624, 629, 604,<br>606, 609, 612, 1109,<br>1112, 1118, 3210                                                                                                                                                                                          |                                                                                                           | ✗             | ✓             | ✓       | ✓        | ✓   | ✓       |
| Microdenier                 | TX59, 3059, 3052                                                                                                                                                                                                                                           | AlphaMops:<br>TX7118M,<br>STX7118M,<br>TX7114M,<br>STX7114M<br><br>BetaMops:<br>TX7070, STX7070           | ✓             | ✓             | ✓       | ✓        | ✓   | ✓       |
| Polyester                   | TX1010, 1012, 1029,<br>1050, 1052, 1060,<br>1069, 1070, 1080,<br>8659, 1003, 1004,<br>1009, 1009B, 1013,<br>1008, 1008B, 2064,<br>2069, 2424, 2452,<br>2409, 2412, 2418, 49,<br>42, 29, 22, 3042, 3049,<br>3215, 3225, 3220,<br>3211, 3212, 3224,<br>8942S | AlphaMops:<br>TX7118, STX7118,<br>TX7114, STX7114<br><br>BetaMops:<br>TX716R, STX716R,<br>TX7072, STX7072 | ✓             | ✓             | ✓       | ✓        | ✓   | ✓       |
| Polyester/Rayon             |                                                                                                                                                                                                                                                            | BetaMops:<br>TX7073, STX7073                                                                              | ✗             | ✓             | ✓       | ✓        | ✓   | ✓       |
| Nylon                       | TX4004, 4009, 4012                                                                                                                                                                                                                                         |                                                                                                           | ✗             | ✓             | ✗       | ✗        | ✗   | ✓       |
| Cotton                      | TX309, TX306, TX304,<br>TX318, TX312, TX312R,<br>TX329, TX3232                                                                                                                                                                                             |                                                                                                           | ✗             | ✓             | ✓       | ✓        | ✓   | ✓       |
| Foam                        | TX704                                                                                                                                                                                                                                                      |                                                                                                           | ✓             | ✓             | ✓       | ✓        | ✓   | ✓       |
| Polypropylene/<br>Cellulose | TX699                                                                                                                                                                                                                                                      |                                                                                                           | ✗             | ✓             | ✓       | ✓        | ✓   | ✓       |

✗ = not compatible    ✓ = compatible



# Pre-Wetted Wipers

| Solution    | Material            | Name         | Size                      | TX#    | Bag Qty | Case Qty              | ISO Class | EU Grade |
|-------------|---------------------|--------------|---------------------------|--------|---------|-----------------------|-----------|----------|
| Non-Sterile |                     |              |                           |        |         |                       |           |          |
| IPA 70%     | Polyester           | Vertex® HS   | 12" x 12" (30 cm x 30 cm) | TX42P  | 50      | 4                     | 3-7       | A-D      |
|             |                     |              | 9" x 9" (23 cm x 23 cm)   | TX49P  | 75      | 4                     | 3-7       | A-D      |
|             |                     | AlphaSat®    | 4" x 4" (10 cm x 10 cm)   | TX1034 | 200     | 4                     | 4-8       | A-D      |
|             |                     |              | 6" x 6" (15 cm x 15 cm)   | TX1036 | 75      | 12                    | 4-8       | A-D      |
|             |                     |              | 9" x 9" (23 cm x 23 cm)   | TX1039 | 50      | 4                     | 4-8       | A-D      |
|             |                     | QuanSat™     | 9" x 9" (23 cm x 23 cm)   | TX1084 | 50      | 12                    | 3-7       | A-D      |
|             | Polyester/Cellulose | TechniCloth® | 6" x 8" (15 cm x 20 cm)   | TX1045 | 100     | 12                    | 5-8       | B-D      |
|             |                     |              | 9" x 11" (23 cm x 28 cm)  | TX1041 | 70      | 12                    | 5-8       | B-D      |
|             |                     |              | 9" x 11" (23 cm x 28 cm)  | TX1065 | 50      | 24                    | 5-8       | B-D      |
|             |                     |              | 7" x 11" (18 cm x 28 cm)  | TX1067 | 200     | 4                     | 5-8       | B-D      |
|             | Polypropylene       | PolySat®     | 7" x 11" (18 cm x 28 cm)  | TX1040 | 200     | 4                     | 5-8       | B-D      |
|             |                     |              | 9" x 11" (23 cm x 28 cm)  | TX1051 | 50      | 24                    | 5-8       | B-D      |
|             |                     |              | 6" x 11" (15 cm x 28 cm)  | TX8723 | 75      | 24                    | 5-8       | B-D      |
|             |                     |              | 6" x 11" (15 cm x 28 cm)  | TX8727 | 75      | 20 & 1 case container | 5-8       | B-D      |
| Ethanol 70% | Polyester/Cellulose | TechniCloth® | 7" x 11" (18 cm x 28 cm)  | TX1068 | 25      | 20                    | 5-8       | A-D      |

|             |                     |              |                           |         |    |    |     |     |
|-------------|---------------------|--------------|---------------------------|---------|----|----|-----|-----|
| Sterile     |                     |              |                           |         |    |    |     |     |
| IPA 70%     | Polyester           | Vertex® HS   | 12" x 12" (30 cm x 30 cm) | TX3042P | 25 | 5  | 3-7 | A-D |
|             |                     |              | 9" x 9" (23 cm x 23 cm)   | TX3049P | 25 | 5  | 3-7 | A-D |
|             |                     | AlphaSat®    | 12" x 12" (30 cm x 30 cm) | TX3252  | 25 | 5  | 4-8 | A-D |
|             |                     | AlphaSat® 10 | 12" x 12" (30 cm x 30 cm) | TX3280  | 50 | 5  | 2-7 | A-D |
|             |                     |              | 9" x 9" (23 cm x 23 cm)   | TX3285  | 20 | 20 | 2-7 | A-D |
|             | Polyester/Cellulose | TechniCloth® | 9" x 11" (23 cm x 28 cm)  | TX3214  | 50 | 24 | 5-8 | B-D |
|             |                     |              | 9" x 11" (23 cm x 28 cm)  | TX3217  | 20 | 24 | 5-8 | B-D |
|             | Polypropylene       | PolySat®     | 9" x 11" (23 cm x 28 cm)  | TX3213  | 50 | 24 | 5-8 | B-D |
|             |                     |              | 9" x 11" (23 cm x 28 cm)  | TX3216  | 20 | 24 | 5-8 | B-D |
| Ethanol 70% | Polyester/Cellulose | TechniCloth® | 7" x 11" (18 cm x 28 cm)  | STX1068 | 50 | 20 | 5-8 | B-D |
|             | Polyester           | Vertex® HS   | 12" x 12" (30 cm x 30 cm) | TX3044P | 25 | 5  | 3-7 | A-D |

# ISO 5 Sample Cleaning Frequency

| Surface             | Each Shift | Daily | Weekly         | Monthly | Quarterly |
|---------------------|------------|-------|----------------|---------|-----------|
| Trash               | ✓          |       |                |         |           |
| Gowning room        | ✓          |       |                |         |           |
| Floors              | ✓          |       |                |         |           |
| Equipment           | ✓          |       |                |         |           |
| Furniture           | ✓          |       |                |         |           |
| Doors               |            | ✓     |                |         |           |
| Windows             |            | ✓     |                |         |           |
| Walls               |            |       | twice weekly ✓ |         |           |
| Ceiling             |            |       |                | ✓       |           |
| Under raised floors |            |       |                |         | ✓         |

Source: IEST-RP-CC-018.4, "Cleanroom Housekeeping: Operating and Monitoring Procedures," p. 13, Table 1.

"A risk assessment should be performed to determine the appropriate frequency for the user. This table is an example of the frequency of cleaning for an average ISO Class 5 cleanroom operation."

## USP<1072> Rotation Recommendations

"Because it is theoretically possible that the selective pressure of the continuous use of a single disinfectant could result in the presence of disinfectant-resistant microorganisms **in a manufacturing area, in some quarters, the rotation of disinfectants has been advocated.**"

"**The daily application of sporicidal agents is not generally favored** because of their tendency to corrode equipment and because of the potential safety issues with chronic operator exposure."

"It is prudent to augment **the daily use of a bactericidal disinfectant with weekly (or monthly) use of a sporicidal agent.**"





Putting the **clean** in **cleanroom** for 50 years.

In 1964 our founder, Edward Paley, created a solution to a problem that no one could see – microcontamination. With the invention of the world's first low-lint wiping cloth, Mr. Paley and Texwipe created an entire industry – contamination control. Today we continue our legacy of clean innovation with products including wipers, swabs, mops, disinfectants, stationery, adhesive mats, sterile products and much more.



Wipers



Swabs



Mops



Disinfectants



Solutions



Stationery



Mats



Sterile



[texwipe.com](http://texwipe.com) | 336.996.7046 | 800.839.9473



|                         |                         |
|-------------------------|-------------------------|
| <b>Texwipe Americas</b> | <b>+1 336 996 7046</b>  |
| <b>Texwipe Europe</b>   | <b>+31 88 1307 410</b>  |
| <b>Texwipe Asia</b>     | <b>+63 049 530 3390</b> |
| <b>Texwipe China</b>    | <b>+65 6468 9433</b>    |
| <b>texwipe.com</b>      |                         |